Chipsmall Limited consists of a professional team with an average of over 10 year of expertise in the distribution of electronic components. Based in Hongkong, we have already established firm and mutual-benefit business relationships with customers from,Europe,America and south Asia,supplying obsolete and hard-to-find components to meet their specific needs.

With the principle of “Quality Parts,Customers Priority,Honest Operation,and Considerate Service”,our business mainly focus on the distribution of electronic components. Line cards we deal with include Microchip,ALPS,ROHM,Xilinx,Pulse,ON,Everlight and Freescale. Main products comprise IC,Modules,Potentiometer,IC Socket,Relay,Connector.Our parts cover such applications as commercial,industrial, and automotives areas.

We are looking forward to setting up business relationship with you and hope to provide you with the best service and solution. Let us make a better world for our industry!

Contact us

Tel: +86-755-8981 8866 Fax: +86-755-8427 6832

Email & Skype: info@chipsmall.com Web: www.chipsmall.com

Address: A1208, Overseas Decoration Building, #122 Zhenhua RD., Futian, Shenzhen, China

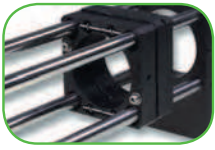
LINOS CATALOG

Edition λ

An Excelitas Technologies Company

Mechanics

High-precision opto-mechanics for reliable integration of optics

Mikrobench
11

Nanobench
79

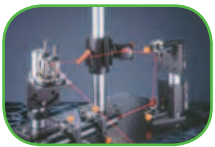
Tube System C
101

Positioning Systems
127


Profile and Rail Systems
167

Mounts, Posts and Holders
189

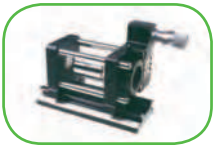Mirror Mounts
223

Q-Sets
237

Optics

Optics with superior performance

Singlets
251

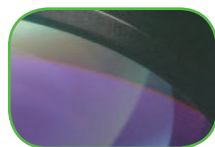
Achromats and Lens Systems
283


Plano Optics
295

Mirrors
331

Thin Film Coatings
359

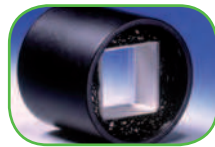Polarization Optics
389

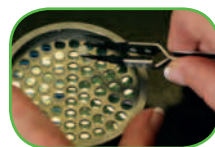
Laser Optics
405

Machine Vision Solutions
423

Micro Imaging Solutions
449

Capabilities
473

Optics Software

Optic Design Software and lens configuration

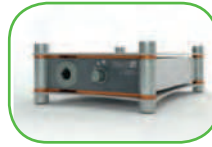
WinLens™ and MachVis
485

Instruments

High-quality laser and light sources for demanding applications

Laser
559

Light Sources
579

Electro Optics

High-performance Faraday isolators, Pockels cells and Modulators

Faraday Isolators
505

Laser Modulators
and Pockels Cells
529

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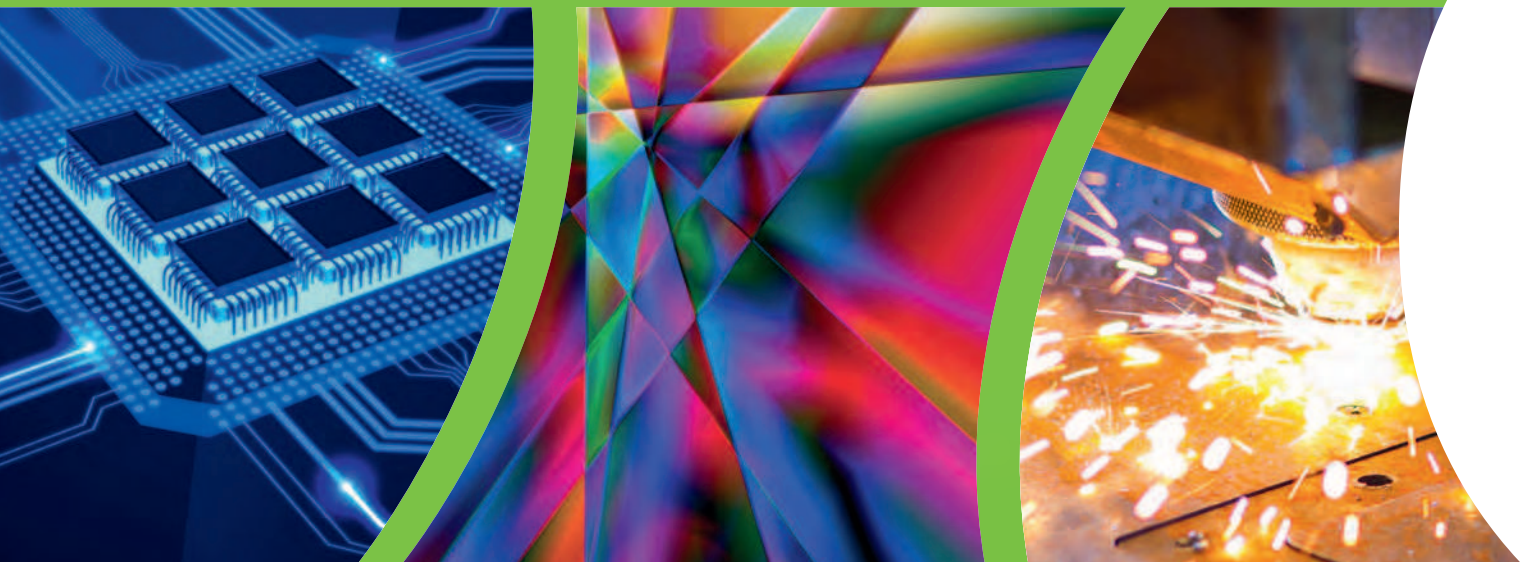
LINOS - Premium brand quality criteria

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Isabel Haackert, Sales Director Catalog & Technical Sales

Dear Readers,

With the new edition λ of the LINOS Catalog, we offer you our extensive range of high-quality standard products. Specifications and useful additional information will help you find solutions for your scientific research requirements, image processing, laser material processing, laser sources and LED illumination and optical measurement technology.

In addition, as a manufacturer of high-end optics we also offer you customized solutions. With our extensive optical fabrication capabilities and design expertise (see page 473), we offer a wide range of optics, from special optics to micro-optics, from aspheres to free form surfaces.

*Use our expertise to meet your requirements! We look forward to receiving your inquiries and being of assistance.
Best regards,*

Isabel Haackert, Sales Director Catalog & Technical Sales

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Excelitas Technologies Corp. is a global technology leader focused on providing innovative, high-performance photonic solutions. Whether you seek off-the-shelf photonic products, or complex custom components, modules and systems, our photonic technologies represent the highest caliber of quality, reliability and performance.

With the acquisition of Qioptiq in 2013, we possess a distinctive brand heritage encompassing pedigrees such as LINOS, Rodenstock, Spindler & Hoyer, Gsanger, Optem, X-Cite, Point Source and OmniCure, amongst others. Our collective knowledge and expertise delivers sophisticated illumination, optical, electro-optical and detection solutions to global customers across the biomedical, scientific research, safety and security, consumer products, semiconductor, industrial manufacturing, and defense and aerospace sectors. Our 5,500 employees operate from 21 state-of-the-art photonics facilities based in 15 different countries across North America, Europe and Asia, and stand committed to fully enabling our customers' innovation and success.

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Visit our online Q-Shop and stay up-to-date! At qioptiq-shop.com you will find up-to-date information about new products, prices and availability. All orders for stock items from Germany that are received working days by 1:00 pm will be sent to you on the same day! You can also request quotes online.

In the Q-Shop you will also find our trade fair and event dates as well as extensive information and numerous publications for download!

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The dialog with our customers is the key to success. Qioptiq employees across three continents are very happy to support you with their expertise and knowledge.

Your Qioptiq contacts

Qioptiq Europe

Phone: +49 (0) 551 6935-0
sales.qioptiq@excelitas.com

Qioptiq North America

Phone: +1 800 775 6786
sales.qioptiq@excelitas.com

Qioptiq Asia Pacific

Phone: +65 6499 7777
sales.qioptiq@excelitas.com

Japan

Opto Science Inc.
Phone: +81 3 3356 1064
info@optoscience.com

Netherlands

VADENO Nederland
Phone: +31 610 414 067
info@vadeno.nl

Mechanics

High-precision opto-mechanics for reliable
integration of optics

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Our quality criteria:

- German manufacturing expertise using only high-quality, extremely durable materials
- Surface finishing with precision anodizing
- Narrow production tolerances
- On-going system expansion with new optical and mechanical components
- The Qioptiq guarantee: Complete compatibility of all Microbench series components - from 1963 to today

LINOS Microbench

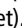
It is our pleasure presenting You the original Microbench, invented by Qioptiq!

This compact 4-rod construction system is in serial production in Göttingen, Germany since 1963. The LINOS Microbench was developed by LINOS predecessor Spindler & Hoyer as an innovative replacement of the triangular rail system. Small in its dimensions, precisely centered, simple to integrate, easy to expand, and very rugged: the LINOS Microbench is ideal for all spatial structures in any compact form.

Thanks to the broad range of optical and mechanical components available, the LINOS Microbench is an integral part of many optomechanical instruments. Take advantage of more than 50 years of experience to ensure the highest precision and stability for your requirements.

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Special Features:

For a rapid setup of your optics, Qioptiq offers a large array of mounted optics developed specially for the Microbench; look for the  symbol in the catalog.

Ideal Areas:

Optics with Ø of 18 to 31.5 mm; illumination and imaging assemblies; optical inspection, measurement and testing equipment; interferometric applications, and more.

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Content

Technical Features	16
Sets	
Microbench Set Mechanics	19
Microbench Set Optics	21
Microbench Basic Set	22
Microbench Set O-UV	24
Laserset	25
Storage Cases	25
Microbench Mechanics Test Kit	26
Mounting Plates	
Mounting Plates 25/30/35/1 "	27
Mounting Plate 25 with Mounting Holes	28
Mounting Plate 25/30/35, Latchable	28
Mounting Plate 25/30, Latchable, 10°	29
Mounting Plate 25 - T10	30
Mounting Plate 25E/30E	30
Mounting Plate M35x0.5, also Latchable	31
Magnetic Cage Assembly Swapper Plates	31
Holder	
Holder 25/30/35/M35x0.5	32
Adjustable Holder 25	32
Rod Holder G/F/90°/V	33
Mounting Bases	34
Rotating Holder, 5 Positions	34
Rods and Cubes	
Rods	35
Cube	35
Beam-Deflector Cube	36
Corner Connectors and Angular Adjustments	
Corner Connector 40/80/40 with Threads	36
Corner Connector T	37
Connector 80 V	37
Swivel Connector	38
Swivel Arm with Scale	38
Mounting Bracket 40	39
Positioners	
XYZ-Fine Adjustment Unit	39
Slip Plate Positioner	40
Centering Mounting Plate/Holder 25	40
X-Y Precision Translation Stage	41
fle.X-Plate	42
Z-Adjustment M	43
Z-Fine Adjustment M	44
Linear Stages for Microbench	45
Z-Translation Stage with Micrometer	46
Z-Translation Stage with Fine-Adjustment Screw	46
Rotary Support	47
Rotary Support 20°	47
Motorized Rotation Mount	48
Counter Support	49
Holder for Micrometer	49
Compression Spring, 60x8x0.7	50

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Content

Fine Adjustment Units and Object Slides

Adjustable Prism Insert 30	50
Adjustable Mirror Insert 30	51
Tilting Platform	51
Prism Table 25, Adjustable	52
Mounting Tool for Beamsplitter Cubes	52
Rotation Adjustment for Cube Inserts	53
Positioning Ring 30	53
Beam Steering Mirror Assembly	54
Optic Support 25	54
Clamp Holder for Prism	55
Prism Support 25	55
Prism Table 25	56
Adjustable Optic Holder 31.5	56

Adapter

Overview Table Adapter	57
Microbench - Nanobench Adapter	58
Adapter 16/25	58
Adapter Collar	59
Cap 30/25	59
Adapter 30/25	59
Adapter 35/30	60
Adapter M35x0.5/30	60
Adapter 35/30-3.5°	60
Adapter 30/C-Mount	61
Adapter 25/C-Mount	61
C-Mount-Adapter-M17	61
Adapter C-Mount/Microbench	62
Adapter 1" to W0.8" or Ø 25 mm	62
Adapter Bench Systems 30 & 1.035"-40 UN	63
Adapter 1.035"-40UN / 30	63
Set of Adapter	64
Adapter 16	64
Adapter 19	64
Adapter 23.2	65
Adapter W0.8"x1/36	65
Adapter W0.8" to 1"	65
Adapter 23.2 to Ø 25	66
Eyepiece Adapter 23.2/30	66
Adapter 25/15	66
FC/FSMA Fiber Adapter	67
FC Fiber Adapter 3.5°	67
Instrument Flange	68
Adapter Plate	68
Set of Retaining Rings	69
Set of Collars	69

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Content

Experiment Set	
Two Beam Interferometer	70
Accessories	
Collimator Holder	71
Cuvette Holder	71
Mounting Plate FLS	72
Closing Caps	72
Focusing Collimators	73
Tube	74
Lens Turret 4x with C-Mount Adapter	74
Microbench Alignment Plate	75
Tools and Screws	
Circlip Kit	75
Special Screwdriver	75
Screwdriver	76
Set of Threaded Pins M2.3x3	76
Set of Threaded Pins M2.3x6	77
Set of Thumbscrews M2.3x9	77
Set of Thumbscrews M2.3x4	77
Set of Screws M2.3x5	77
Set of Hexnuts M2.3	78
Polyamide Plugs	78

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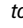
Technical Features

Microbench is a compact component system for numerous areas of application – from lighting and imaging structures to optical experimental setups, measurement / inspection equipment and interferometric applications, to name just a few. The basic principle lies in the combination of standardized mounting plates for integrating optics of all types, connected by stainless steel rods. Both linear and spatial setups can be implemented quickly and easily. All components of the Tube Mounting System C with outer diameter of 30 respectively 35 mm can easily be integrated in the appropriate Microbench mounting plates. Via adapters the LINOS Microbench is also compatible with our Nanobench system.


Qioptiq offers mounted optics developed specially for the Microbench with diameters from 18 mm to 31.5 mm. The special  symbol shows you at a glance which optics are available for your Microbench system.

- Construction system for
 - illumination or imaging systems
 - optical testing, metrology, or inspection systems
 - interferometric systems
 - Compact, three-dimensional setups
- Rigid 4-rod structures
- Optional 2-rod structures for more rapid setup
- Easy to handle
- High mechanical stability

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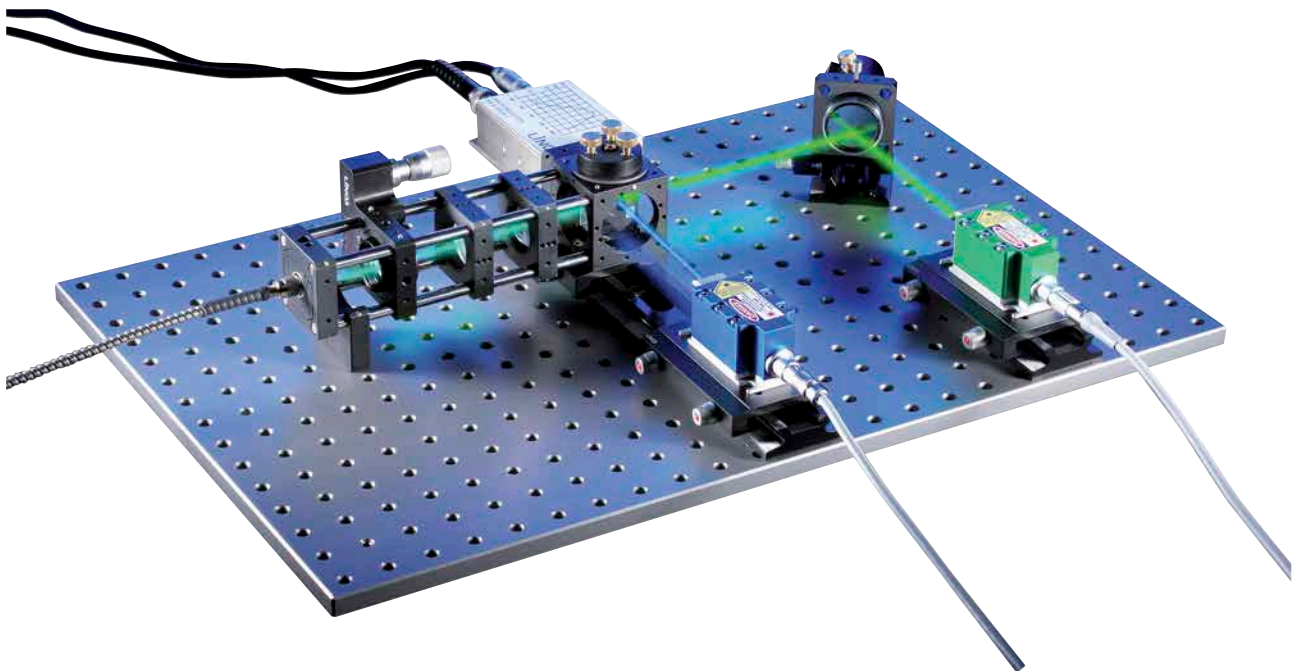Please find further information about application examples, guidelines, frequently asked questions and more in our Brochure "Microbench, Bench and Rail systems - Introduction, solutions and applications". You will find the brochure for download at each Microbench product under "Docs + Drawings".

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Mounting plates

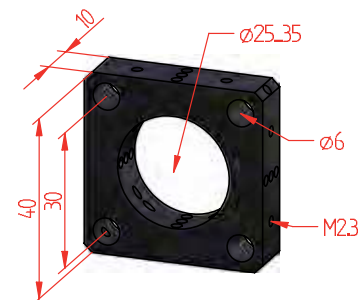
- Lead-free aluminum alloy
- Matt black anodized surfaces
- Inner diameter: 25, 30 or 35 mm with H7 tolerances
- Four bore holes for mounting the connecting rods. The tolerances for the rod bore holes on the mounting plates, together with the tolerances for the rods, guarantee a system with virtually zero play.

Rods


- Hardened stainless steel with polished surfaces
- Lapped to fit
- Outer diameter: 6 mm
- Lengths: 20 mm to 450 mm
- Compatible adjustment rings as holders or stops; e.g., for Z-adjustment of individual components

Cubes

- Multifunctional elements for 3D structures
- Ideal for beam deflectors
- Inner diameter of bore holes: 25 or 30 mm
- Two of the cube sides have bore holes for rods, four sides have threads for mounting plates with tapped holes

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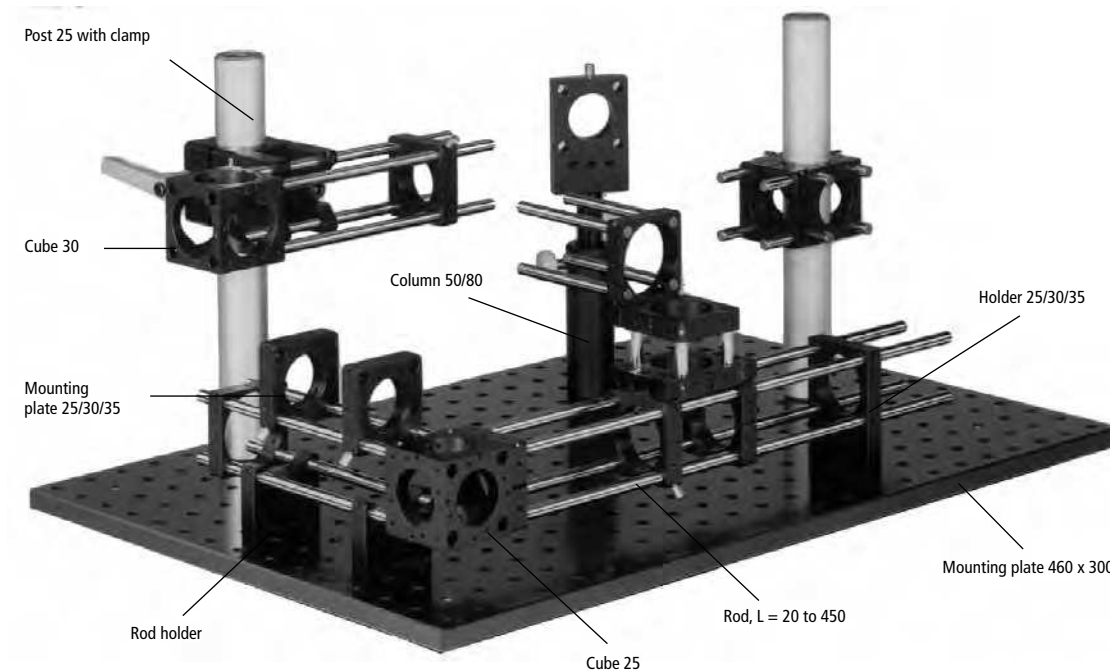- Includes all most important parts for an easy start with Microbench system
- Complete with rails and carriers of the Rail System FLS 40
- Contains tools and installation material

Microbench Set Mechanics

Part No.	Quantity	Description
G06 1355 000	1	Rail FLS 40-250
G06 1374 000	1	Carrier FLR 40-40
G06 1372 000	1	Carrier FLR 40-20
G06 1371 000	2	Carrier FLR 40-10
G06 1309 000	1	Mounting Plate 220x40 mm
G06 5061 000	2	Holder 25
G06 1225 000	2	Holder 30
G06 5084 000	2	Holder 35
G06 1010 000	2	Mounting Plate 25
G06 1042 000	3	Mounting Plate 30
G06 1047 000	2	Mounting Plate 35
G06 1020 000	1	Mounting Plate 25, latchable
G06 1021 000	1	Mounting Plate 30, latchable
G06 1041 000	2	Mounting Plate with mounting holes
G06 1025 000	2	Centering Mounting Plate 25
G06 1080 000	1	Cube 25
G06 1081 000	1	Cube 30
G06 1228 000	1	Rod Holder
G06 1229 000	1	Rod Holder G
G06 1227 000	2	Rod Holder F
G06 1110 000	4	Corner Connector 40 mm
G06 1161 000	1	Counter Support
G06 1160 000	1	Holder for micrometer G 10
G06 1162 000	1	Micrometer G 10
G06 1159 000	2	Compression Spring
G06 5058 000	1	Rotary Support
G06 3730 000	1	Adjustable Mirror Insert 30
G06 3731 000	1	Adjustable Prism Insert 30
G06 5002 000	1	Optic Support 25
G06 1661 000	1	Adapter 1" to W 0.8 and ø 25 mm
G06 1601 000	1	Cap 30/25
G06 1621 000	1	Adapter W0.8x1/36"
G06 1636 000	1	Adapter 23.2
G06 1602 000	1	Adapter 30/25
G06 1658 000	1	Adapter 30/C-Mount
G06 1637 000	1	Adapter 35/30
G06 1650 000	1	Iris Diaphragm C 18 with tube 25
G06 1220 000	1	Set of 10 collars
G06 5092 000	1	Positioning Ring 30
G06 1207 000	4	Rod, 20 mm
G06 1208 000	4	Rod, 30 mm
G06 1209 000	4	Rod, 75 mm
G06 1210 000	4	Rod, 150 mm
G06 1205 000	4	Rod, 200 mm

Microbench Set Mechanics

Item Title	Part No.
Microbench Set Mechanics	G060004000

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- Fast beginning of experiments ensured by wide range of mounted optics

Microbench Set Optics

Part No.	Quantity	Description
G06 3020 000	1	Biconvex lens 20/21.4 mounted
G06 3024 000	1	Biconvex lens 50/21.4, mounted
G06 3026 000	1	Biconvex lens 80/21.4, mounted
G06 3028 000	1	Biconvex lens 150/21.4, mounted
G06 3852 000	1	Biconvex lens 50/24, mounted
G06 3854 000	1	Biconvex lens 100/24, mounted
G06 3014 000	1	Biconvex lens 70/30, in mount 31.5
G06 3042 000	1	Plano-convex lens 30/17, mounted
G06 3043 000	1	Plano-convex lens 40/21.4, mounted
G06 3044 000	1	Plano-convex lens 50/21.4, mounted
G06 3045 000	1	Plano-convex lens 60/21.4, mounted
G06 3046 000	1	Plano-convex lens 80/21.4, mounted
G06 3047 000	1	Plano-convex lens 100/21.4, mounted
G06 3048 000	1	Plano-convex lens 150/21.4, mounted
G06 3049 000	1	Plano-convex lens 200/21.4, mounted
G06 3821 000	1	Plano-convex lens 50/24, mounted
G06 3822 000	1	Plano-convex lens 80/24, mounted
G06 3823 000	1	Plano-convex lens 100/24, mounted
G06 3826 000	1	Plano-convex lens 300/24, mounted
G06 3052 000	1	Plano-convex lens 100/30, in mount 31.5
G06 3057 000	1	Biconcave lens -20/21.4, mounted
G06 3060 000	1	Biconcave lens -40/21.4, mounted
G06 3062 000	1	Biconcave lens -100/21.4, mounted
G06 3892 000	1	Biconcave lens -50/24, mounted
G06 3070 000	1	Plano-concave lens -40/21.4, mounted
G06 3871 000	1	Plano-concave lens -50/24, mounted
G06 3072 000	1	Plano-concave lens -100/21.4, mounted
G06 3076 000	1	Concave-convex lens 50/21.4, mounted
G06 3127 000	1	Achromat 40/17, mounted
G06 3141 000	1	Achromat 50/21.4, mounted
G06 3212 000	1	Achromat 60/24, in mount 25.4
G06 3143 000	1	Achromat 80/21.4, mounted
G06 3144 000	1	Achromat 100/21.4, mounted
G06 3215 000	1	Achromat 120/24, in mount 25.4
G06 3204 000	1	Achromat 160/30, in mount 31.5
G06 3148 000	1	Achromat 200/21.4, mounted
G06 3191 000	1	Achromat -50/17, mounted
G06 3010 000	1	Condenser combination 16/21.4
G06 3099 000	1	Asph. condenser lens 18/21.4, mounted
G06 3098 000	1	Asph. condenser lens 27/30 in mount
G06 3450 000	1	Colored glass filter BG 38, mounted
G06 3451 000	1	Colored glass filter BG 18, mounted
G06 3082 000	1	Colored glass filter VG 9, mounted
G06 3081 000	1	Colored glass filter RG 630, mounted
G06 3080 000	1	Heat absorbing filter KG 1, mounted
G06 3410 000	2	Polarization filter
G06 3411 000	1	Quarter-waveplate
G06 3619 000	1	Front surface concave mirror 14/30 mounted
G06 3628 000	1	Front surface plane mirror ø 30 mounted
G34 0060 000	1	Front surface plane mirror 20 x 30 mm
G34 4143 000	1	Beamsplitter plate 20 x 30 mm
G39 0137 000	1	Plane plate 20 x 30 x 2.5 mm
G33 9913 000	1	90°-Prism 20 x 20 mm
G33 5520 000	1	Beam splitter cube 20 x 20 x 20 mm
G03 8771 000	1	Microscope objective 10x/0.25

Microbench Set Optics

Item Title	Part No.
Microbench Set Optics	G062004000

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Microbench Basic Set

The case contains all components of the following sets:

Microbench Set Mechanics G060004000
Microbench Set Optics G062004000

- Types of systems that may be assembled:
 - Microscopes
 - Off-axis viewers
 - Telescopic systems
 - Measuring telescopes
 - Collimators
 - Polarizers
 - Projection systems
 - Graticule projectors
 - Beamsteering systems
 - Beamsplitter systems
 - Interferometers
 - Illumination systems
- Included: Software Package WinLens™ (see chapter Optics Software for further information)

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Please find further information about application examples, guidelines, frequently asked questions and more in our Brochure "Microbench, Bench and Rail systems - Introduction, solutions and applications".

You will find the brochure for download at each Microbench product under "Docs + Drawings".

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
Microbench Set O-UV

Item Title	Part No.
Microbench Set O-UV	G062020000

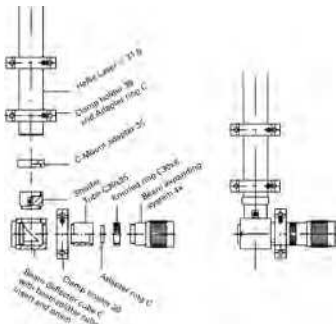Mounting example:
Horizontal angle assembly with beam expander system

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Storage Cases

All storage cases complete with inserts, but without components.

- Suitable for careful and organized storage of Microbench components and components with similar dimensions
- Layout of the compartments as shown for the sets with components

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